



① PREPARE


SYRINGE WITH 0.5M  $\text{NiCl}_2$  SOLUTION AND WITH DISTILLED WATER  
5% RUBEANIC ACID (YOU WILL NEED THAT IN ABOUT TWO DAYS)  
VASELINE SYRINGE

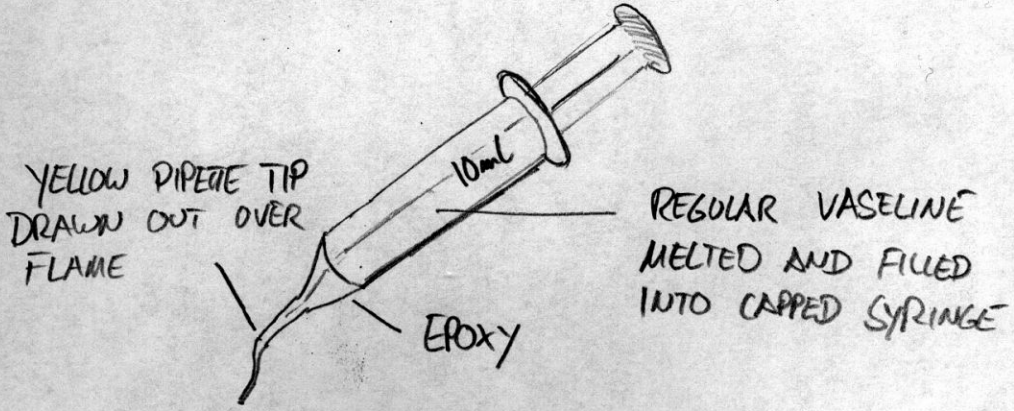



② LAY BOTTOM PLATE UNDER CLEANLY CUT NERVE ENDING

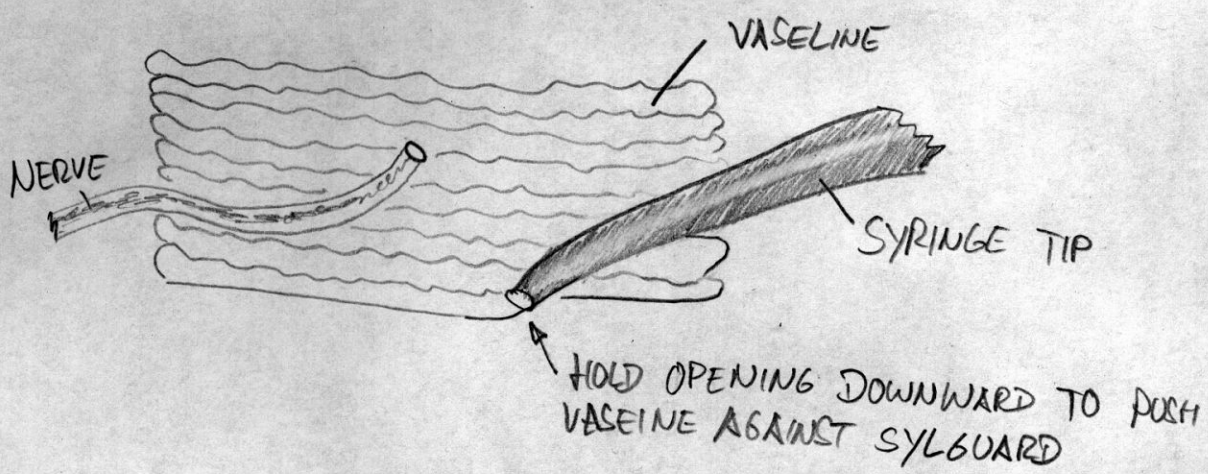

③ BUILD UP WELL, RING BY RING

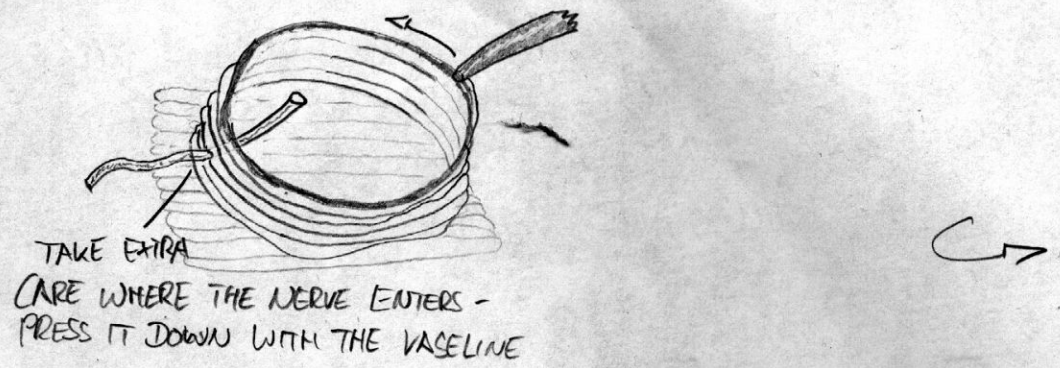


- ④ WHEN THE WELL BREAKS THE SALINE LEVEL, TEST IF IT IS WATER TIGHT

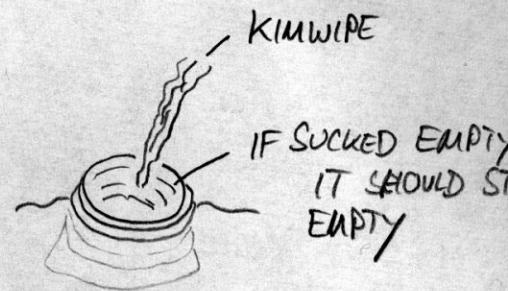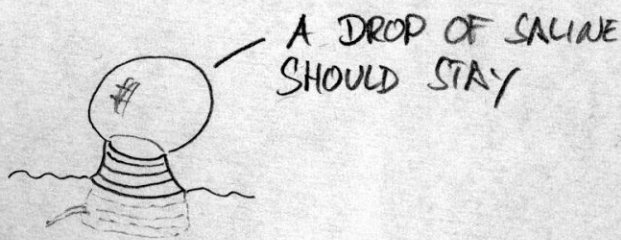

- ⑤ SUCK SALINE FROM WELL AND FILL WITH A DROP OF DISTILLED WATER

- ⑥ AFTER 1 MINUTE SUCK OUT DISTILLED WATER AND FILL WITH  $\text{NiCl}_2$ .

- ⑦ COVER WELL

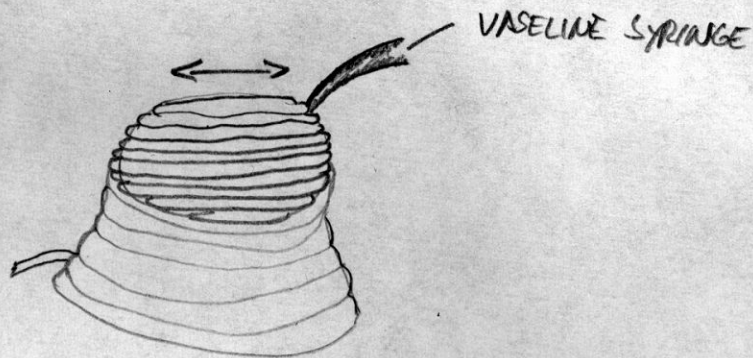

- ⑧ PUT PREP DISH IN A LARGER CONTAINER WITH DAMP TISSUE AND PLACE IN COLDROOM ( $4^\circ\text{C}$ ) FOR  $\sim 36$  HOURS

- ⑨ CAREFULLY REMOVE COVER WITH FORCEPS, SUCK OUT  $\text{NiCl}_2$ , LOWER SALINE LEVEL IN PREP DISH. DROP 5% RUBENIC ACID ONTO PREP. AFTER 5 MINUTES YOU SHOULD SEE BLACK NERVES. REMOVE REST OF WELL AND WASH WITH SALINE.